

# I N S T R U C T I O N S



## Maintenance of Scienceware® Eye Wash Bottles

Catalog No. F24850-0000, F24851-0000

1. Wash the Eye Wash Bottle and all accessory parts: cap, tubes, etc. in mild detergent solution.
2. Rinse well to remove all traces of detergent.
3. Under clean conditions, fill Eye Wash Bottle completely.

*(We recommend Scienceware® sterile eye wash solution refill # F24879-0032. Please note that once the refill bottle has been opened, the sterility of the solution cannot be assured).*

4. Replace top, close cap and return to emergency station.
5. If an eye wash bottle is used, empty it completely and proceed through steps 1-4 before replacing in station.
6. After using an eye wash bottle, the injured individual should see a physician at once for examination and treatment.

To ensure cleanliness, it is advisable to establish a regular eye wash bottle cleaning and eye wash solution replacement procedure. See the refilling guidelines below.



## Guidelines for Refilling Scienceware® Eye Wash Bottles

The below recommendations for solution replacement are based on information gathered from sources outside Bel-Art Products to help assist our customers. Bel-Art Products does not guarantee the accuracy of the information gathered.

Rinse solution and frequency of change:

- *Water only – change once a week.*
- *Water and Preservative (varies by manufacturer) - change every two months.*
- *Pre-mixed Eye Wash Solution - change every 4 to 6 months.*
- *Unopened, Sterile Eye Wash Solution (F24879-0032) - until expiration date on bottle.*

Bel-Art Products Sterile Eye Wash solution (F24879-0032) has a shelf life of up to 3 years - refer to expiration date on bottle. However, once the bottle has been opened, the product is no longer sterile. The Sterile Eye Wash Solution is meant for use with Scienceware® Eye Wash Stations (F24878-1032 and F24878-2032).

If using the Sterile Eye Wash Solution (F24879-0032) to refill Scienceware® Emergency Eye Wash Bottles (F24850-0000 and F24851-0000), the solution can no longer be considered sterile once transferred into those bottles. Replace transferred solution every 4 to 6 months.

**BEL-ART PRODUCTS**

PEQUANNOCK, NJ 07440-1992 USA  
TEL: 1-800-4BEL-ART • FAX: 973 694-7199 • www.belart.com